PrimerCube®

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OPERATING INSTRUCTIONS

Please read the operating instructions before using this device.

Only use the device as described in these operating instructions.

Keep these instructions for future reference.

PrimerCube®

PACKAGE CONTENTS:

1x Primer Cube rotary magnetic oscillator
1x USB-C charging cable
1x 5V 2A USB charging adapter
1X operating instructions

PRIMERCUBE FUNCTION:

The rotary magnetic oscillator is functionally adjustable with the high accuracy of the selected speed. Choice of speed in the interface: 900–3500 Rpm/min
Note: Rpm – Revolutions per minute

The device has a built-in button for setting the application time; the application time can be selected in three time intervals.

Warning: The PrimerCube device is sensitive to shocks. It has a built-in motor with a rotating pan equipped with permanent neodymium magnets. This pan rotates at high speed. Make sure to avoid any blows and shocks!!!

The purpose of the device: laboratory use, physical experiments with a magnetic rotating field.

BATTERY CHARGING PRECAUTIONS

The device batteries must be fully charged before using the PrimerCube for the first time. When charging the battery for the first time, charge the PrimerCube for at least 8–10 hours. Subsequent full charging of the battery will take approximately 3-5 hours. The type of battery built in PrimerCube: 8,8 A / h Li-ION. This type of battery has no memory function. Consequently the battery can be charged at any capacity phase, it is not necessary to discharge the battery completely in order to be charged.

Discharging the battery completely shortens the battery life!

ONLY USE THE ORIGINAL CHARGER SUPPLIED WITH PRIMERCUBE!
It is recommended to use PrimerCube with

a connected charger. This reduces the internal batteries wear thus extending the life of PrimerCube significantly.

We only recommend using PrimerCube powered by batteries for a short time (3–5 hours)

The charging proceeds as follows: after connecting the USB charging cable and the adapter to the power supply, connect PrimerCube to the charger. The charging process is indicated by a red control light located next to the USB connector. The first charging should take at least 8-10 hours. The display shows the battery status and the charging process status. The battery charging process is fully automatic, and the charging process ends automatically when the battery is fully charged.

We recommend using PrimerCube during charging when possible. It can also be used temporarily with an internal battery.

DISPLAY FUNCTIONS AND DEVICE OPERATION

1: Display activation

Connect the original charger to the USB power connector on the PrimerCube and charge it for at least 8–10 hours.

Hold down the START = / Zzz / button on the upper left corner of the display for 5–7 seconds. When pressed properly, the red LED will start flashing – wait until the display lights up.


After this, the display leaves the STANDBY mode and is displayed with fully functional buttons.

2: Warning

Before using the PrimerCube / battery for the first time or in an event the battery has not been used for a long time, it is recommended to charge it to the full (8-10 hours of charging without operation). Only use chargers / adapters and batteries approved by the manufacturer. During the charging process, the display shows the battery symbol and its status; the status of the charging process is displayed next to this marking. The red LED next to the USB connector lights up.

Before using for the first time, set the current time in the upper right corner by touching and then using the numeric keypad. Thanks to the correctly set time, the device can correctly apply the timer function (T1, T2, T3).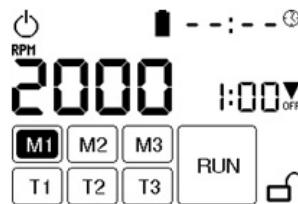


3: Application time setting

Before the device is used set the application time.

Touch the M1 key on the display. Use this button to set application time No.1. Then press the button showing the application time (on the right in the middle of the display time arrow OFF) and then the display with numerical options will appear:

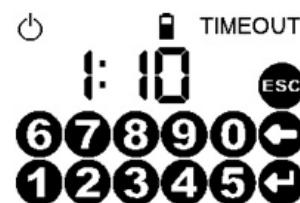After setting the required application time press the ENTER – button This saves the settings.

Choose the speed after this step.

4: Setting M1, M2, M3 – memory (speed, application time)


The RPM and application time below are set in the factory settings of the device memory:

M1: 1,000 RPM 2 hours
M2: 2,000 RPM 2 hours
M3: 3,350 RPM 1 hour

These factory settings can be changed according to individual needs.

5: RPM speed setting

Touch the display to set the desired RPM speed. After pressing the numeric RPM value, the following numeric keypad will appear on the display: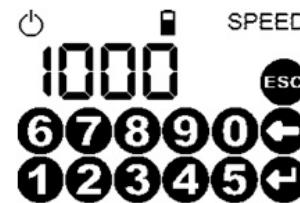


Use this numeric keypad to set the desired RPM speed.

Press ENTER to confirm and save the required settings.

6: Change of the application time default setting

The application time is shown on the right in the middle section of the display. Touch this displayed application time to set the desired application time. After pressing this time value, the following numeric keypad will be displayed: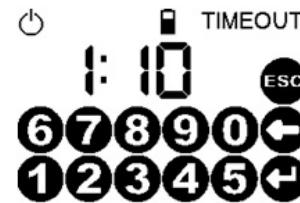

Use this numeric keypad to set the desired application time.

Press ENTER to confirm and save the required settings

Setting and saving individual programs is possible at positions M1, M2, M3; you can use them to choose your own required values of the application time and RPM speed values.

7: Setting and using the T1, T2 and T3 timer


A program with an individual timer can be set in positions T1, T2, T3. Thus, the PrimerCube will operate in a preset time interface.

Note: Example: PrimerCube can be pre-programmed at the position T1, T2 or T3 to be activated and then deactivated in the time interval from 03:00 to 07:00 a.m. The device is switched on at 3.00 a.m. and switched off at 7.00 a.m. automatically.

Press the T1 key:
The option to set two time values appears in the upper right corner of the display. These two time values show the beginning and end of the duration of the set program:

These two time values can be set and saved by touching and then using the numeric keypad: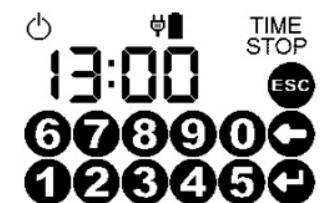

The timer can be activated with the EN button. After pressing the EN button, you can see that the T1 position is displayed in a solid dark rectangle. This means that the set timer is **active** and the PrimerCube is **switched on and off** at the required time.

The deactivation of this function – timer – can be set by pressing the DIS button. After deactivation, the T1 value is displayed without a border in a smaller black cube.

After the time interval of 5 minutes elapses, during which the PrimerCube has not been active, the device switches to the STANDBY mode, i.e. the sleep mode. In this off mode, the internal battery can be maintained fully charged for about 1 month. Preset programs are saved in the device's memory because the EEPROM memory does not depend on the battery charging activity.

If the battery level drops below 5 %, the PrimerCube will automatically turn off and OFF will appear on the display. In this case, connect the PrimerCube to the power supply source using a charging cable / adapter and charge for at least 8–10 hours.

A button with the „lock“ symbol is located in the lower right corner of the display. Use this button to LOCK. When this function is active, the touch screen closes securely and no settings can be changed on it; neither incidentally nor intentionally.

8: On/Off

After activating the desired program by pressing the RUN button, the in-built rotary motor spins. The active operation of the device motor is indicated on the display by the M letter in circle. At the same time, during the operation of the device, the in-built red LED above the display lights up. Thanks to this, it is easy to check the device activity even in entire darkness.

LEDs control the luminous intensity according to the ambient light value. Thus, when in complete darkness, the luminosity of LEDs dims to a minimum within a few minutes after start-up so that the user is not disturbed, for example during sleep.

Note Example: After starting the M1 program, the display will show the RPM speed value and the remaining application time until the end of the program is displayed to the right of the RPM speed. After this application time elapses the PrimerCube is switched off. The picture shows the remaining application time: 58 minutes. The LOCK function is not active. The lock is open.

If the user wants to stop the program before the remaining application time has elapsed, press the STOP button on the display.

TECHNICAL PARAMETERS

- Battery: capacity 8.8Ah, Li-ION.
- 2-level protection against overcharging and discharging.
- Short circuit protection and over-voltage protection.
- RPM: 900–3,500
- The electronics are secured against overheating.
- Regulation of light intensity according to the surroundings.
- Connector: USB-C max 12V / 5A.
- USB-C standards: PD3.0 / PD2.0 / QC3.0 / QC2.0 / FCP / PE2.0 / PE1.1 / SFCP.
- Adapter: QC3.0, 12V / 1.5A.
- The device can be used as a PowerBank in an emergency.
- Motor: max 3,500rpm, 30W.

TROUBLESHOOTING:

PrimerCube does not work, the display does not respond – connect the PrimerCube to the original charger (regardless of the battery status on the display) and charge at least 3 hours. Then use a paper clip – there is a small hole on the PrimerCube next to the charging connector – the reset button. Press the reset button for 3 seconds using a paper clip with caution. Then your PrimerCube gets to the factory settings and everything should work smoothly again. Then charge your PrimerCube batteries with the original charger to 100%

For other problems, contact our technical support: info@primercube.eu

FIRE SAFETY

If your device is extremely hot or smells, do not attempt to remove the battery. If you detect smoke coming from your device, clear the area around it. Sprinkle the fire with sand or use an approved fire extinguisher, or cover it with a fireproof container. Never use water to extinguish, as it may react with lithium and cause a fire. If you can move the device safely, place it in a fireproof or well-ventilated area and wait until it cools completely. Therefore, in case of any doubts, turn off the device, place it in a fireproof container and get it repaired in an authorized repair shop.

WARNINGS AND PRECAUTIONS

Use the device correctly and safely - therefore,

read these warnings and precautions carefully and follow them while using the device. Unless stated otherwise, the term "device" means the device and its accessories.

When storing, carrying and using the device, keep it dry and protect it from collision with other objects. Do not disassemble the device. In an event of a malfunction, contact an authorized service repair shop for assistance or repairs. Neither any organization or an individual is allowed to change the mechanical, safety and performance parameters of the device without permission. When using the device, comply with all applicable laws and regulations and respect the rights of others.

Requirements for the environment, in which the device is used. Keep the device away from heat and fire sources such as heaters or candles. Keep the device away from electronic devices that produce strong magnetic or electric fields, such as microwave ovens or refrigerators. Place the device on a stable surface. Keep the device in a cool and well-ventilated indoor environment. Never expose the device to direct sunlight. During charging, the device should be in an ambient temperature of 0°C to 35°C. When using a battery-powered device, the ambient temperature should range within 0°C and 35°C. Do not place any objects (such as a candle or water container) on the device. If an object or liquid enters the device, stop using the device immediately, turn it off, disconnect all connected cables, and then contact an authorized repair shop.

During a thunderstorm, turn off the device and then disconnect all connected cables from the device to protect it from a damage caused by the lightning.

SAFETY DURING TRAVEL

Local laws and regulations must be complied with when using the device. In addition, if you use the device while driving, follow these instructions: Focus on driving. Your primary responsibility is to drive safely. Radio frequency (RF) signals can adversely affect electronic systems in motor vehicles. Do not place the device on an air bag or in an area where the air bag is deployed in a motor vehicle. Failure to do so could result in a device damage due

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to excessive force when inflating the air bag. Never use the device while aboard an aircraft. Switch off your device before boarding an aircraft. The use of wireless devices in an aircraft may be dangerous to the operation of the aircraft and disrupt the wireless telephone network. In addition, this might be considered illegal.

Never use the device in cases when the use of wireless devices is prohibited or may cause interference or danger. If you use an electrical medical device, contact the manufacturer of your electrical device for information on restrictions on the device use. Never bring the device to operating rooms, intensive care units or cardiovascular diseases units.

FLAMMABLE AND EXPLOSIVE AREAS

Never use the device at places where flammables or explosives are stored, such as gas stations, oil tanks or chemical plants. Otherwise, there is a risk of explosion or fire. In addition, follow the instructions in the warnings or symbols. Do not store or carry the device in the same box with flammable liquids, gases, or explosives.

Accessory requirements: Only use accessories supplied together with the product or authorized by the device manufacturer. Otherwise, the performance of the device and the device warranty may be affected, or an accident may happen. Do not use the AC adapter if its cord is damaged. Otherwise, there may be a risk of electric shock or fire. Make sure that the power adapter meets the parameters listed on the device nameplate. Make sure that the power adapter meets the requirements of Article 2.5 of the IEC60950-1 / EN60950-1 guideline and that it has been tested and approved in accordance with state or local standards. For devices that can be plugged into an outlet, the outlet must be located near the device and must be easily accessible. Unplug the charger from the electrical plug and the device when not in use. If the power cord is damaged (for example, the damaged insulation or the cord is broken), or the plug is loose, stop using the cord immediately. Failure to do so could result in electric shock, short circuit of the charger or fire. Use the AC power source specified in the charger specifications. Improper power supply

voltage may cause fire or charger failure.

SAFETY OF CHILDREN

Keep the device and its accessories away from children. Otherwise, the device and its accessories could be accidentally damaged, there is a risk of swallowing small parts of the device and suffocation or other dangerous situations.

MAINTENANCE

If you do not use the device for a long time, turn it off and then disconnect all cables connected to it. If an unexpected accident occurs, such as smoke escaping from the device or if the device makes an unusual noise or smells, stop using the device immediately, turn it off, disconnect all connected cables, and then contact an authorized repair shop. If the screen of the device is damaged by a heavy object impact, neither touch it nor attempt to repair the component. In this case, stop using the device immediately and then contact an authorized repair shop. Do not step on, pull on or twist the cord. Failure to comply with this instruction may damage the cable and cause the device malfunction. Before cleaning, stop using the device, turn it off and then disconnect all cables connected to it. Use a clean, soft, dry cloth to clean the surface of the machine. Do not use any cleaning agents or sprays to clean the surface of the device.

WARRANTY CONDITIONS

The company grants the warranty for the period of 24 months from the date of purchase of the product subject to the compliance with instructions for installation, operation and maintenance of the product.

The warranty does not cover any defects caused by unprofessional intervention in the use or neglecting care and maintenance of the device and defects caused by improper device handling contrary to the instructions for use.

Internal battery life is 6 months from purchase/ from commissioning.

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